




止痛药在NanoChrom BP-5MS上的分离与检测

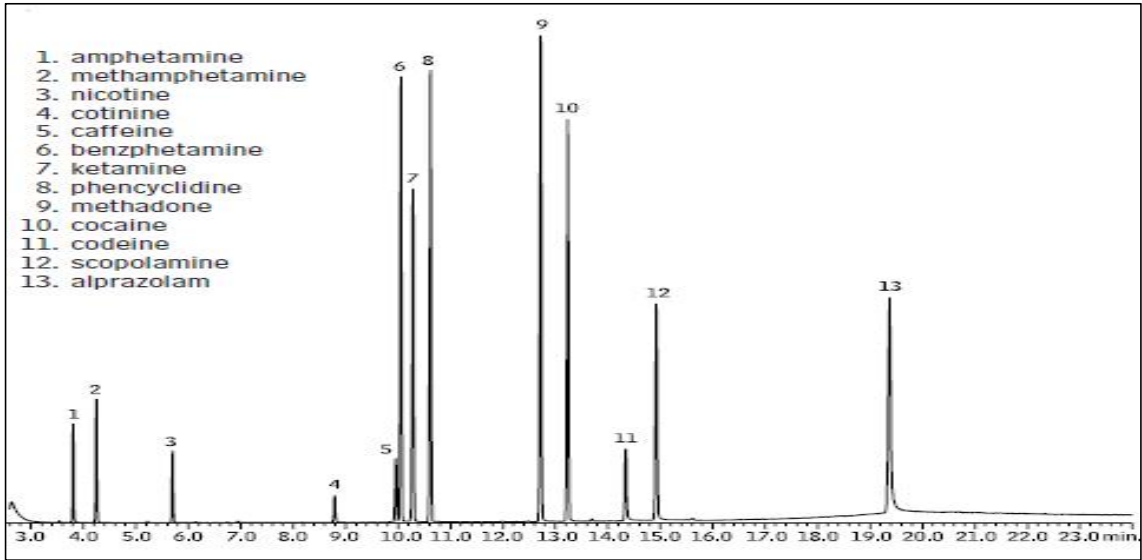

Column: NanoChrom BP-5MS
 Dimension: 30m×0.25mm×0.25µm
 Carrier: He, flow 1.0 mL/min
 Oven: 100 °C to 220 °C to 320 °C (hold 3 min) at 10 °C/min
 Inject: 0.2 µL
 Detector: MSD
 Peaks:

1. amphetamine (安非他明)
2. methamphetamine (甲基苯丙胺)
3. nicotine (尼古丁)
4. cotinine (可替宁)
5. caffeine (咖啡因)
6. benzphetamine (苯丙胺)
7. ketamine (氯胺酮)
8. phencyclidine (苯环定)
9. methadone (美沙酮)
10. cocaine (可卡因)
11. codeine (可待因)
12. scopolamine (东莨菪碱)
13. alprazolam (阿普唑仑)